



## КАПИЛЛЯРНЫЕ КОЛОННЫ DN-RAN

**Архангельск** (8182)63-90-72  
**Астана** (7172)727-132  
**Астрахань** (8512)99-46-04  
**Барнаул** (3852)73-04-60  
**Белгород** (4722)40-23-64  
**Брянск** (4832)59-03-52  
**Владивосток** (423)249-28-31  
**Волгоград** (844)278-03-48  
**Вологда** (8172)26-41-59  
**Воронеж** (473)204-51-73  
**Екатеринбург** (343)384-55-89  
**Иваново** (4932)77-34-06  
**Ижевск** (3412)26-03-58  
**Казань** (843)206-01-48

**Калининград** (4012)72-03-81  
**Калуга** (4842)92-23-67  
**Кемерово** (3842)65-04-62  
**Киров** (8332)68-02-04  
**Краснодар** (861)203-40-90  
**Красноярск** (391)204-63-61  
**Курск** (4712)77-13-04  
**Липецк** (4742)52-20-81  
**Магнитогорск** (3519)55-03-13  
**Москва** (495)268-04-70  
**Мурманск** (8152)59-64-93  
**Набережные Челны** (8552)20-53-41  
**Нижний Новгород** (831)429-08-12  
**Новокузнецк** (3843)20-46-81

**Новосибирск** (383)227-86-73  
**Омск** (3812)21-46-40  
**Орел** (4862)44-53-42  
**Оренбург** (3532)37-68-04  
**Пенза** (8412)22-31-16  
**Пермь** (342)205-81-47  
**Ростов-на-Дону** (863)308-18-15  
**Рязань** (4912)46-61-64  
**Самара** (846)206-03-16  
**Санкт-Петербург** (812)309-46-40  
**Саратов** (845)249-38-78  
**Севастополь** (8692)22-31-93  
**Симферополь** (3652)67-13-56  
**Смоленск** (4812)29-41-54

**Сочи** (862)225-72-31  
**Ставрополь** (8652)20-65-13  
**Сургут** (3462)77-98-35  
**Тверь** (4822)63-31-35  
**Томск** (3822)98-41-53  
**Тула** (4872)74-02-29  
**Тюмень** (3452)66-21-18  
**Ульяновск** (8422)24-23-59  
**Уфа** (347)229-48-12  
**Хабаровск** (4212)92-98-04  
**Челябинск** (351)202-03-61  
**Череповец** (8202)49-02-64  
**Ярославль** (4852)69-52-93

**DN-PAH**  
Dedicated Columns

**DN-PAH 15m**

ID	Film	Max Temp	Code	Chroma
0.25mm	0.25µm	350°C	9414.117 303	067

**DN-PAH 25m**

ID	Film	Max Temp	Code	Chroma
0.25mm	0.25µm	350°C	9414.117 304	


**DN-PAH** Technical Specifications

Every Column Individually Tested  
 Test Certified and Grob Mixture included in each Column  
 Instruction Manual included in each Column

DANI DN-PAH Capillary Column

Bonded and cross-linked  
 Inertness  
 Low bleeding  
 Good thermal stability


Similar to NO EQUIVALENT

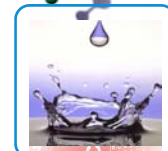

# DN-PAH

## Chromatograms

### Polyaromatic Hydrocarbons (PAHs)

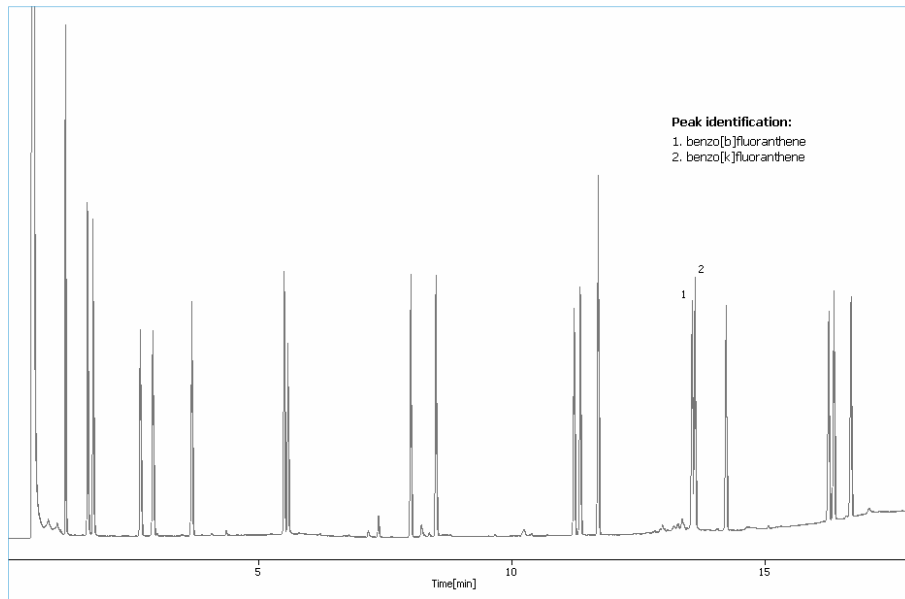

DN-PAH 15m 0.25mm 0.25µm

#### Chromatographic Conditions

<b>Oven</b>	140°C - 1.0 min - 10°C/min 350°C - 1.0 min
<b>Injection</b>	PTV - 40°C - 999°C/min - 400°C - 5 min
<b>Carrier Gas</b>	Hydrogen - 1.0 ml/min
<b>Injection volume</b>	0.5 µL
<b>Detector</b>	FID - 400°C

067



# DN-PAH FAST

Dedicated Columns

## DN-PAH FAST 15m

ID	Film	Max Temp	Code	Chroma
0.10mm	0.10µm	350°C	9414.117 302	066

## DN-PAH FAST Technical Specifications

Every Column Individually Tested  
 Test Certified and Grob Mixture included in each Column  
 Instruction Manual included in each Column

DANI DN-PAH FAST Capillary Column

Bonded and cross-linked  
 Inertness  
 Low bleeding  
 Good thermal stability


Similar to NO EQUIVALENT



# DN-PAH FAST

## Chromatograms

Fast GC application: Polyaromatic Hydrocarbons (PAHs)


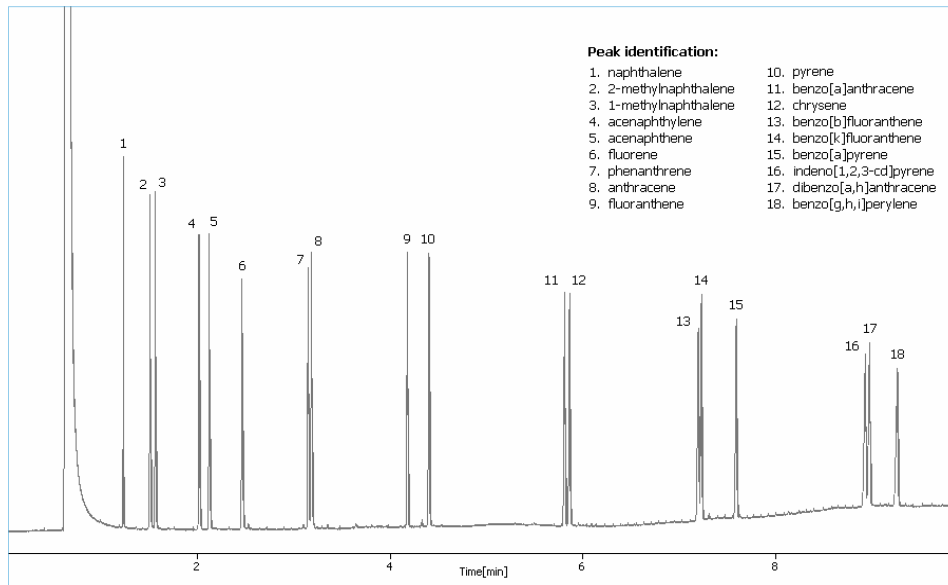


DN-PAH FAST 15m 0.10mm 0.10µm

**Chromatographic Conditions**

<b>Oven</b>	140°C - 0.5 min - 30°C/min 220°C - 15°C/min - 300°C
<b>Injection</b>	PTV - 80°C - 999°C/min - 400°C
<b>Carrier Gas</b>	Hydrogen - 0.5 ml/min
<b>Injection volume</b>	0.5 µL
<b>Detector</b>	FID - 400°C

066

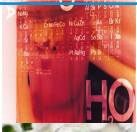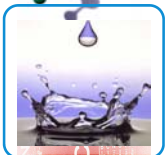

**По вопросам продажи и поддержки обращайтесь:**

**Архангельск** (8182)63-90-72  
**Астана** (7172)727-132  
**Астрахань** (8512)99-46-04  
**Барнаул** (3852)73-04-60  
**Белгород** (4722)40-23-64  
**Брянск** (4832)59-03-52  
**Владивосток** (423)249-28-31  
**Волгоград** (844)278-03-48  
**Вологда** (8172)26-41-59  
**Воронеж** (473)204-51-73  
**Екатеринбург** (343)384-55-89  
**Иваново** (4932)77-34-06  
**Ижевск** (3412)26-03-58  
**Казань** (843)206-01-48

**Калининград** (4012)72-03-81  
**Калуга** (4842)92-23-67  
**Кемерово** (3842)65-04-62  
**Киров** (8332)68-02-04  
**Краснодар** (861)203-40-90  
**Красноярск** (391)204-63-61  
**Курск** (4712)77-13-04  
**Липецк** (4742)52-20-81  
**Магнитогорск** (3519)55-03-13  
**Москва** (495)268-04-70  
**Мурманск** (8152)59-64-93  
**Набережные Челны** (8552)20-53-41  
**Нижний Новгород** (831)429-08-12  
**Новокузнецк** (3843)20-46-81

**Новосибирск** (383)227-86-73  
**Омск** (3812)21-46-40  
**Орел** (4862)44-53-42  
**Оренбург** (3532)37-68-04  
**Пенза** (8412)22-31-16  
**Пермь** (342)205-81-47  
**Ростов-на-Дону** (863)308-18-15  
**Рязань** (4912)46-61-64  
**Самара** (846)206-03-16  
**Санкт-Петербург** (812)309-46-40  
**Саратов** (845)249-38-78  
**Севастополь** (8692)22-31-93  
**Симферополь** (3652)67-13-56  
**Смоленск** (4812)29-41-54

**Сочи** (862)225-72-31  
**Ставрополь** (8652)20-65-13  
**Сургут** (3462)77-98-35  
**Тверь** (4822)63-31-35  
**Томск** (3822)98-41-53  
**Тула** (4872)74-02-29  
**Тюмень** (3452)66-21-18  
**Ульяновск** (8422)24-23-59  
**Уфа** (347)229-48-12  
**Хабаровск** (4212)92-98-04  
**Челябинск** (351)202-03-61  
**Череповец** (8202)49-02-64  
**Ярославль** (4852)69-52-93

**Единый адрес для всех регионов: [drs@nt-rt.ru](mailto:drs@nt-rt.ru) || [www.danimaster.nt-rt.ru](http://www.danimaster.nt-rt.ru)**

